



KERN & Sohn GmbH

Ziegelei 1
72336 Balingen-Frommern
Germany

+0049-[0]7433-9933-0
FAX +0049-[0]7433-9933-149
info@kern-sohn.com

Operating manual KERN Interface Adapter with Cable for Bluetooth

KERN YKUP-06

Typw TYKUP-06-A
Version 1.0
2021-09
GB



You will find the current version of these instructions also online under:
<https://www.kern-sohn.com/shop/de/DOWNLOADS/>
Under the column Operating instructions

TYKUP-06-A-BA-e-2110_Bluetooth



KERN Interface Adapter with Cable

Version 1.0 2021-09

Installation Instructions for Bluetooth

Contents

1	General hints	2
1.1	Installation.....	2

1 General hints

Cable length: 0.15 m

i	<ul style="list-style-type: none">• Weighing data can be transferred via Bluetooth.• Only KERN KUP-adapters may be connected to the 15-pol-sub-D-connection of the balance!
----------	---

1.1 Installation

- Switch off appliance
- Plug in KUP-adapter (Bluetooth) at the 15-pol-Sub-D-connection of the appliance
- Switch on appliance

The KUP adapter is optionally available with **Bluetooth Low Energy** (BLE) and will then be visible to Bluetooth Master devices using its serial number.

For access, please use an appropriate software application / app which supports Bluetooth Low Energy (BLE). Applications exclusively using Bluetooth Classic (BTC) will not work.

The following profile must be set:

Service UUID
0000fff0-0000-1000-8000-00805f9b34fb

Read characteristic UUID
0000fff1-0000-1000-8000-00805f9b34fb

Write characteristic UUID
0000fff2-0000-1000-8000-00805f9b34fb